

FLUORESCENT PROTEIN IMAGING

TYPHOON USE FOR DIGE IMAGING

(SE 600 gels)

1. Clean Platen with water
 - a. Use water ONLY.
 - b. Do not open gels, wash and dry.
2. Place the Black Gripper labeled “Front” at the bottom of the platen (the side nearest you.)
3. Place the Black Gripper labeled “Back” at the top of the platen (the side farthest from you).
4. Place the Black Gripper labeled “Center” in the middle of the platen. The two ends should fit snugly into the grooves centered on both the “Front” and “Back” grippers.
5. Place the small black grippers on the left and right sides of the platen. The groove on each gripper should face inwards and two grippers should be placed on each side.
 - a. This allows four gels to be scanned at one time.
 - b. Holding the gel in its normal orientation, slide one side of the gel into the groove of a small black gripper, and place the opposite side on the metal tabs protruding from “Center” gripper.
 - c. It’s best to practice putting the gel on the scanner the same way every time.

